

02-JAN-1997

LABEL (1)

MICROTABS(TM)

used to preserve the integrity of laboratory milk samples

35,000 TABLETS

READ LABEL / PACK INSERT BEFORE USING

KEEP OUT OF THE REACH OF CHILDREN

Store in a dry place, between 15°C and 30°C

(TM) 'MICROTABS' is a trade mark of D & F Control Systems, Inc.

GUARANTEE: Contains 54% 2-bromo-2-nitropropane-1,3 diol
(equivalent to 10mg active per tablet)

REGISTRATION NO. 21508 PEST CONTROL PRODUCTS ACT

D & F Control Systems, Inc.
3401 Crow Canyon Road, Suite 110, San Ramon,
California, 94583, U.S.A.
Tel.: 510 830 0429
Fax.: 510 830 0379

AGENT:

P.A.T.L.Q., Dairy Herd Analysis Service
MacDonald College, Poultry Building
Ste.-Anne de Bellevue, Quebec, Canada H9X 1C0
Telephone: 514-398-7880

DANGER

POISON

EYE IRRITANT

CORROSIVE

FIRST AID: In the case of personal contact with skin or eyes
give prolonged irrigation with water. In the case of ingestion
wash out the mouth thoroughly with water and give water to drink.
In all cases obtain medical attention.

PRECAUTIONARY STATEMENTS: Severe eye or skin irritant. Avoid
ingestion, inhalation and contact with skin, eyes and clothing.

Where there is risk of contact, wear protective clothing.

Store away from: (a) Oxidising materials. (b) Heat and sources of ignition. (c) Food and food containers. Avoid contact with metals.

TOXICOLOGICAL INFORMATION: The tablets and concentrated solutions are skin irritants.

LABEL (2)

MICROTABS(TM)

used to preserve the integrity of laboratory milk samples

COMMERCIAL

READ THE LABEL / PACK INSERT BEFORE USING

DANGER

POISON

EYE IRRITANT

CORROSIVE

GUARANTEE: Contains 54% 2-bromo-2-nitropropane-1,3 diol

REGISTRATION NO. 21508 PEST CONTROL PRODUCTS ACT

CONTAINS: 800 TABLETS IN EACH DISPENSER

PATENTED DISPENSER - FOR EXCLUSIVE USE WITH DFC PRODUCTS

KEEP OUT OF THE REACH OF CHILDREN

Store in a dry place between 15°C and 30°C

D & F Control Systems, Inc.
3401 Crow Canyon Road, Suite 110, San Ramon,
California, 94583, U.S.A.
Tel.: 510 830 0429
Fax.: 510 830 0379

((PAMPHLET))

MICROTABS(TM)

COMMERCIAL

READ THE LABEL BEFORE USING

REGISTRATION NO. 21508 PEST CONTROL PRODUCTS ACT

READ THE LABEL BEFORE USING

DANGER
EYE IRRITANT

POISON
CORROSIVE

D & F Control Systems, Inc.
3401 Crow Canyon Road, Suite 110, San Ramon,
California, 94583, U.S.A.
Tel.: 510 830 0429
Fax.: 510 830 0379

Guarantee: Each tablet contains 54% 2-bromo-2-nitropropane-1,3 diol (equivalent to 10mg active per tablet).

Use Level: One tablet per 50 mls (2 fl. ounces) of milk sample.

Solubility: At room temperature the active is fully dispersed within 15 minutes if use directions on dispenser are followed.

Safety: The active is not sensitizing to healthy human skin and is not irritant at normal usage levels. For further information on the toxicity, consult D & F Control Systems, Inc.

Storage: Tablets should be stored in a dry place between 15°C and 30°C. Keep away from food and food containers, sunlight and fluorescent lights. Keep out of the reach of children.

Biological Activity: Active against bacteria but not yeast or mould. Milk samples must be adequately refrigerated during transportation and storage at the laboratory in order to prevent growth of yeast or mould. Samples held between 5° C (40° F) and 10° C (50°

F) during transportation and storage should be tested within three days. Samples held under 5°C (40° F) during transportation and storage should be tested within five days.

Colour

Indicator: MICROTABS contain a colour indicator that allows detection of spoiled samples. A normal colour of blue-grey indicates that the sample is satisfactory for testing. A yellow milk sample indicates spoilage and should not be tested. A blue-grey sample having a yellow surface layer (fat layer) indicates spoilage from yeast or mould and should not be tested.

Laboratory

Instrument

Calibration: The use of MICROTABS results in minimal interference when using normal laboratory techniques. To obtain maximum instrument accuracy when testing milk samples, this product may be included in calibration and control milks at normal laboratory usage levels.

SPILLAGE: Immediately sweep up and remove, then clean the area with detergent and water.

WASTE DISPOSAL: Burn or bury in compliance with local regulations.

NOTICE TO USER: This control product is to be used only in accordance with the directions on this label. It is an offence under the Pest Control Products Act to use a control product under unsafe conditions.

NOTICE TO BUYER: Seller's guarantee shall be limited to the terms set out on the label and subject thereto, the buyer assumes the risk to persons or property arising from the use or handling of this product and accepts the product on that condition.

™ 'MICROTABS' is a trademark of D & F Control Systems, Inc.

KEEP OUT OF THE REACH OF CHILDREN

For further information contact:

Mr. Brian Corrigan, Lab Manager, P.A.T.L.Q.
Dairy Herd Analysis Service,
MacDonald College, Poultry Building,
Ste.-Anne de Bellevue Quebec, Canada H9X 1C0

Telephone: 514-398-7880

This label transcript service is offered by the Pest Management Regulatory Agency to provide efficient searching for label information. This service and this information do not replace the official hard-copy label. The PMRA does not provide any guarantee or assurance that the information obtained through this service is accurate, current or correct, and is therefore not liable for any loss resulting, directly or indirectly, from reliance upon this service.

+))